# Autoloop PRO

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**Flame time**  
*Precise settings of the required flame sterilization time*

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**Cooling period**  
*The cooling time can be set for the inoculation loop material used - exact to the second*

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**Sterilization monitor**  
*An intelligent sensor monitors the safe sterilization. Three setting options / off switch*

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### Easy to use

*All parameters are easy to set; quick start button for immediate start*

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### Mount safely

*Coordinated mounts and reliable optical detectors for inoculation loop holders*

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### Optimal sterilization

*Sterilization monitor for safe sterilization of inoculation loops*

\* Fußpedal optional / Foot pedal optional

# Autoloop<sup>PRO</sup>

## The fully automatic carousel for flame sterilizing inoculation loops.

The inoculation loop carousel **Autoloop<sup>PRO</sup>**, which is suitable for up to 4 inoculation loops, was developed to comply with current requirements in microbiological, cellbiological and biotechnological laboratories.

### Modern design

The stable housing enables comfortable and easy access to inoculation loops. Removal positions on both sides make the carousel equally suitable for right and left handers.

### More comfort

Keep all functions in view with the fully graphic display. The graphic menu control makes it easy to select all functions quickly. Flaming and cooling time can be adjusted to the second. During the flame sterilizing, the display shows the remaining flaming time. The **Autoloop<sup>PRO</sup>** rotates and controls flaming automatically.

No unintentional use of hot inoculation loops: When flaming is completed, the display shows the remaining cooling time and the removal positions of cool inoculation loops. All the suspended inoculation loops can be flamed as needed in one full rotation of the carousel.

For additional flame sterilization (e.g. of bottle necks or test tubes) a foot pedal\* can be activated during the automatic operation. For this purpose the **Autoloop<sup>PRO</sup>** turns the inoculation loop out of the flame position and enables additional flame sterilization.

### More safety

The nearly perpendicular position of the inoculation loop holder protects even the sleeve nut of the holder from possible contaminations. Additionally, an intelligent sensor of the **Autoloop<sup>PRO</sup>** monitors safe sterilization.

### Saves energy, time and costs

Precise flame sterilization time extends the life of the inoculation loops and reduces gas consumption. Continuous working during the flaming and the cooling phases of the inoculation loops saves a great deal of time and makes it possible to work efficiently.

### Robust

Fabricated entirely of stainless steel, anodized aluminium and a display, protected by heat resistant glass, the **Autoloop<sup>PRO</sup>** can withstand extreme laboratory conditions.

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### Start options

Start immediately or sterilize all inoculation loops in a full rotation

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### Everything visible at a glance

Removal positions of cool inoculation loops or the remaining cooling time, residual heat display and sterilization time

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### NEW:

**Additional flame sterilization** of bottle necks or test tubes during the automatic operation by activating a foot pedal.

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## The range:

**Autoloop<sup>PRO</sup>**

includes 4 inoculation loop holders  
4 inoculation loops  
Pedestal  
Foot pedal connector  
Instruction manual  
2-year warranty  
Laboratory gas burner **not**  
included

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**Spray protector for Autoloop<sup>PRO</sup>**  
for more safety when handling with  
pathogenic material

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**Spare inoculation loops**  
made of special stainless steel,  
wire Ø 0.6 mm

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CE: EN 61326-1, EN 61010  
EWG Richtlinien: 2004/108/EG, 2006/95/EG

## Technical data:

Technology: Microprocessor  
Display: fully graphic dot-matrix  
Rotational direction: left / right, depends on removed holder  
Inoculation loop holder: 1 - 4, with optical holder detection  
Time of flame sterilizing: 1 - 15 sec  
Time for cool-down period: 1 - 45 sec  
Additional flame sterilization: by foot pedal (optional)  
Safety features: temperature control  
flame activation warning  
indicator of remaining cooling time  
tilt sensor, residual heat display

Adjustable  
sterilization monitor: 1 - Off  
2 - low intensity  
3 - middle intensity  
4 - high intensity

Stand by: automatic unit switch off: 1 h  
(or with stand-by time of the burner)

Operating voltage: via gas burner (foot pedal socket)  
Casing: stainless steel & aluminum,  
UV and solvent resistant

Measurements: Pedestal (Art.-No. 8.000.400):  
135 x 180 mm (w x d)  
Pedestal (Art.-No. 6.000.400):  
135 x 135 mm (w x d)

Height Autoloop<sup>PRO</sup>: 305 mm  
Weight: Art.-No. 8.000.400: 1200 g  
Art.-No. 6.000.400: 1100 g

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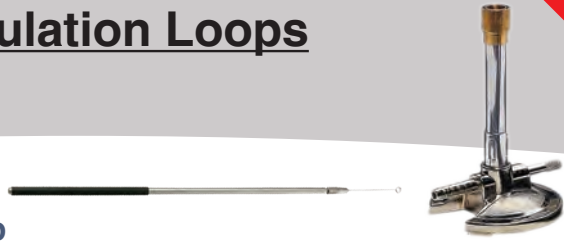
**SAVE UP TO  
70% TIME**

## Speed-up Flame Sterilizing Inoculation Loops and Cut Down Your Costs!

### Case 1: Bunsen Burner & 1 Inoculation Loop

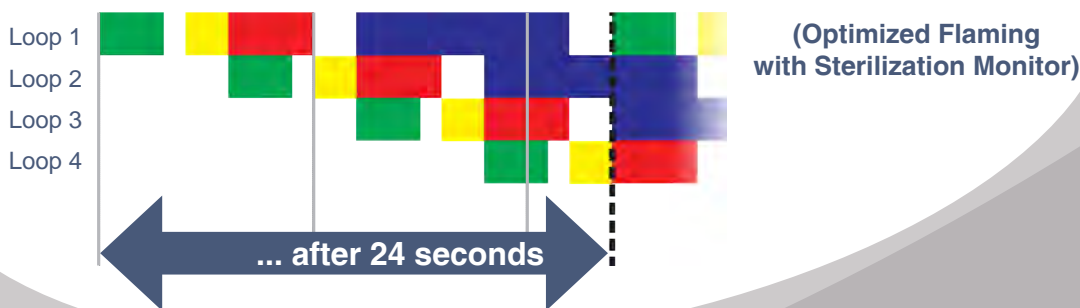
- Working with inoculation loop
- Flaming inoculation loop
- Time to change loops
- Cooling down period

### Case 2: Fuego SCS & 4 Inoculation Loops placed on a Tray

### Fall 3: Fuego SCS with Automated Carousel Autoloop Pro & 4 Inoculation Loops

**Working with the Autoloop Pro (Case 3):**  
 Save 70% working time compared to Case 1  
 Save 40% working time compared to Case 2

**Compared to Case 2 the Autoloop Pro has paid off after approx. 50 working hours!\***

\* based on an hourly rate of 30 Euro